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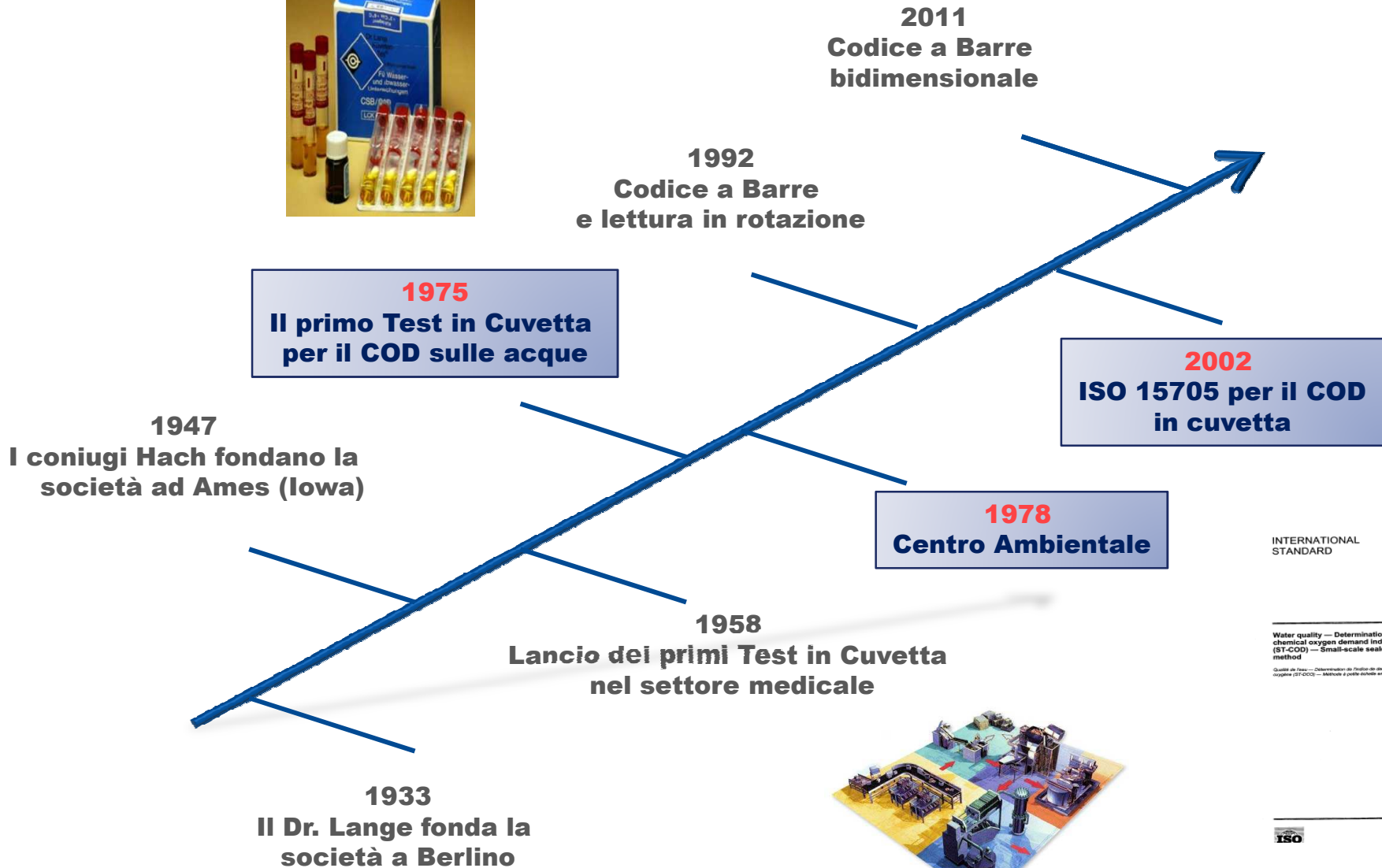
Hach Lange Srl
Stefano Malusardi
Lab Manager

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
INTERNATIONAL STANDARD
ISO 15705
First edition
2002-11-15

Water quality — Determination of the chemical oxygen demand index (ST-COD) — Small-scale sealed-tube method

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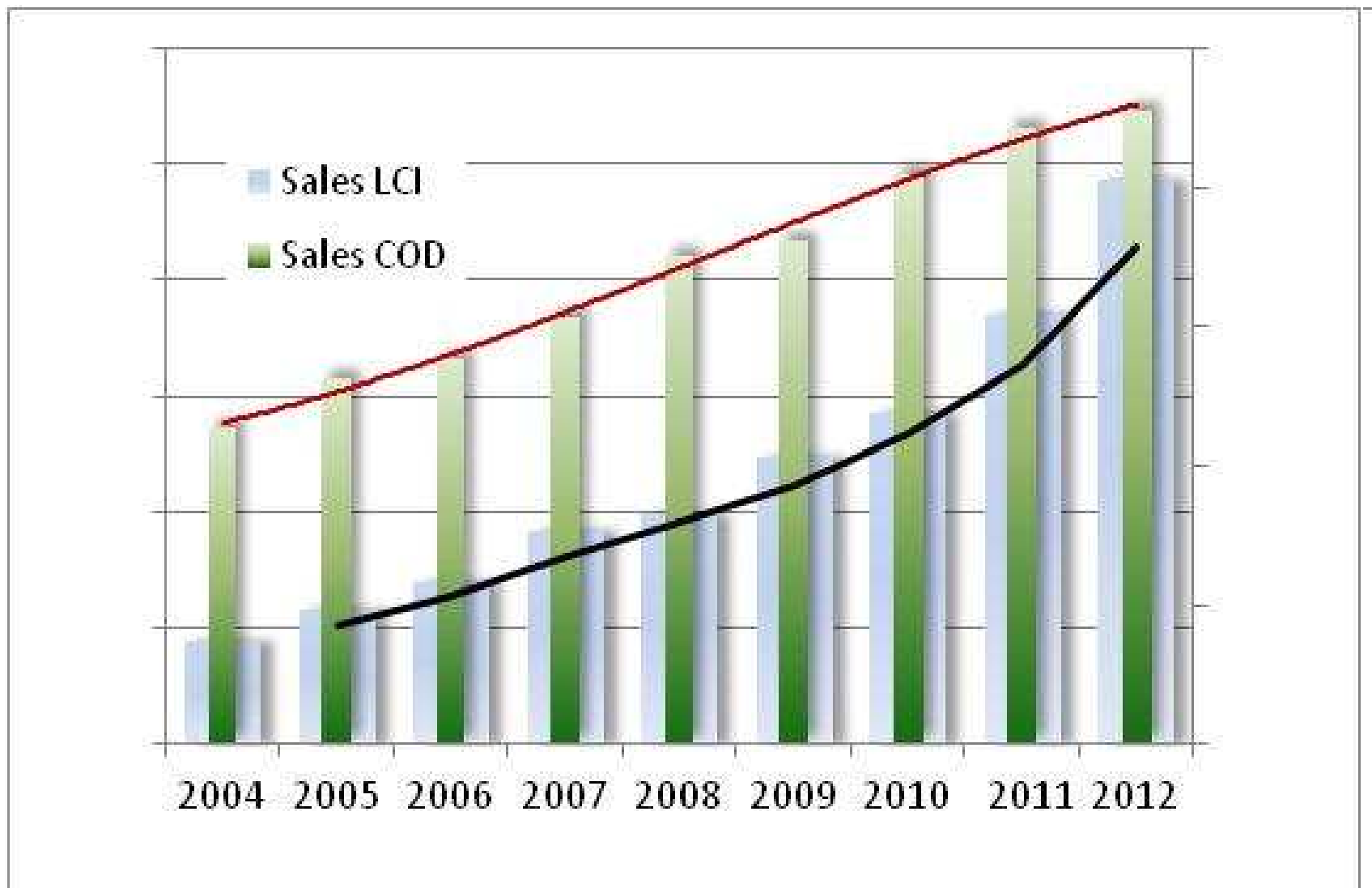Certificate of Conformity

We hereby declare that the
LANGE Cuvette Test LCI 500
 is completely conform to
ISO 15705
 „Determination of the chemical oxygen demand index (ST-COD)-
 Small-scale sealed tube method“

Quality certificate
 Technical data for Validation
 of LI500 (0-150 mg/l COD in
 conformity with ISO 15705)

Required
 Formulae
 one: K₁C
 one: K₂C
 one: K₃C
 one: K₄C

Sample V
 ml: 20 ml

Digestion
 Temperature
 Time: 2 h

Probiostat
 Wavelength

HACH LAM
 L500 Appl
 Düsseldorf

Dr. Ralf

HACH LANGE
 UNITED FOR WATER QUALITY

Quality certificate
 Technical data for cuvette test LCI500

Sensitivity	0.0032 Abs./mg/l
Ordinate intercept	0.003 Abs.
Residual standard deviation	0.0026 Abs.
Method variation coefficient	0.96 %
Method standard deviation	0.8 mg/l
Confidence interval: (95%)	± 1.9 mg/l

The technical data for cuvette test LCI500 were determined in conformity with ISO 8466-1 and DIN 38402 A51 „Calibration of analysis methods“.

The series of the smallest and largest calibration standards exhibit normal distribution and are outlier- and trend-free.

The calibration gives a linear function.

Technical data in conformity with DIN 38465

Detection limit	2.3 mg/l
Quantitation limit	6.9 mg/l

The detection and the quantitation limits were determined in conformity with DIN 38465.

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ISPRA

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- Collaborazioni con Commissioni Europee:
 - **Environment (ENV)**
 - **Agriculture and Rural Development (AGRI)**
 - **Joint Research Centre (JRC)**
 - **Maritime Affairs and Fisheries (MARE)**
- Iniziative al Parlamento Europeo:
 - **Environment, Public Health & Food Safety**
 - **Internal Market & Consumer Protection**
 - **Regional Development**
 - **Agriculture & Rural Development**

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